

Premium-Level Micro/Semi-Micro Balances

A&D Borealis

BA-T/BA Series



Sophisticated Weighing at a Touch

AND
A&D Company, Ltd.


Discover Precision

www.aandd.jp

Supreme Precision, Optimizing Your Weigh

Come experience premium weighing solutions with outstanding ease of use, realized by A&D Borealis BA-T/BA series. In addition to promising fully reliable results, it provides exceptional efficiency and time savings to improve your daily tasks and routines while meeting various laboratory requirements. An ideal choice if you seek topnotch professional quality and need nothing but the best!

BA-T series

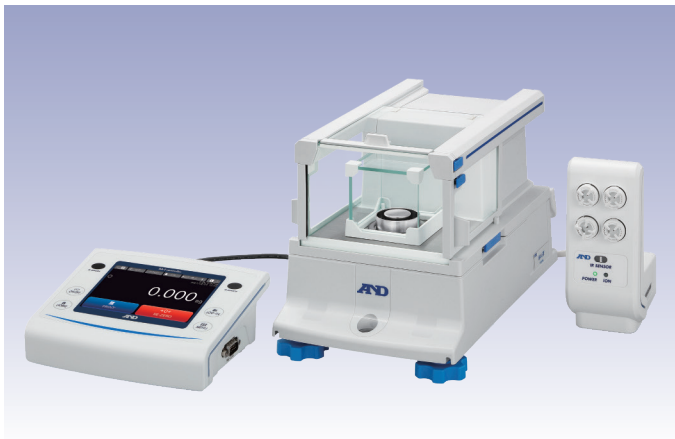

0.001 mg models (micro balances)



0.01 mg models (semi-micro balances)

With a 5-inch wide color touch screen, the BA-T series brings maximum operability, enhanced functionality and a broader range of solutions. All models come standard with an external ionizer (static eliminator).

BA series



0.001 mg models (micro balances)



0.01 mg models (semi-micro balances)

The BA series has a standard reverse-backlit LCD, and is suitable for users who only require core functionality. The 0.001 mg models are equipped with an external ionizer as standard.

✓ Operating a balance should never be a source of distress, even when weighing micrograms

Automatic opening/closing of the breeze break doors

The side doors of the breeze break can be opened/closed using the non-contact IR sensors on the display unit (or optional foot switches) for smooth, efficient access to the weighing chamber. For the actuator, A&D adopt a pump and air cylinder mechanism,♦ which is much more quiet and durable than conventional motor actuators.

♦ Patent pending

Breeze break unit

The top door can be oriented so that the handle is positioned to the rear instead of the front, making it easier for you to see inside.

Inner breeze break for 0.001 mg models♦

The open side can be easily turned around.

♦ Patent pending

Weighing unit

IR sensors

For the BA-T series, the IR sensors can also be used for contactless operation of the RE-ZERO or PRINT command.



Automatic doors

You can set the opening distance to full, half, or your preferred distance.



Touch screen and user navigation (for the BA-T series)

The touch screen enables intuitive operations while making it easy to enter numbers, change settings, etc. You can operate it even if you are wearing thick gloves as it responds to pressure (i.e. resistive touch screen).

Display unit

Four frequently used keys are provided as physical keys for quick access.

✓ Innovative design that makes cleaning the balance simple and effortless

Easily separable glass panes

The each glass pane of the breeze break can be easily separated for thorough cleaning/disinfection thanks to the unique clip system.♦

♦ Patent pending



No grooves for the sliding doors♦

The doors suspend from the ceiling instead of resting on grooves so that contaminants do not accumulate during use.

♦ Patent pending

Detachable breeze break unit

The breeze break unit itself can also be easily removed from the weighing unit by pulling out the two side fasteners. ♦ This allows for fast, simple cleaning as well as replacement in case of failure.


♦ Patent pending



✓ How utmost weighing performance, compliance and security can be assured at all times

On-screen leveling assist (for the BA-T series)

The touch screen provides a visual instruction on which leveling foot (feet) to rotate in which direction(s) depending on the position of the air bubble on the level indicator.

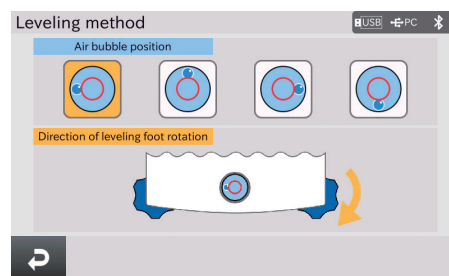

AD-Just – Automatic self sensitivity adjustment

The balance can be set to calibrate and adjust its sensitivity automatically using its internal weight according to the set execution condition (i.e. temperature change, preset time or interval) when there is nothing on its weighing pan. An indicator (message for the BA-T series) blinks to give notice before the adjustment starts.

Internal repeatability test

It is also possible to have the balance test and calculate repeatability (standard deviation) using its internal weight to quickly assess the performance under a given environment.

External ionizer for instant static removal

With the provided ionizer,*1 you can easily ensure that your sample (and container) is completely free from destabilizing static electricity. The ionizer adopts the DC method, which does not require fans to deliver ions, and therefore causes no breeze (except for minimal ionic wind). This allows you to remove static without disturbing even extremely fine powders.

The ionizer draws power via the balance. For seamless weighing workflows, it can be installed next to the balance and activated for a set duration by hovering your hand over the embedded IR sensor right before a sample is placed inside the breeze break.

*1 Optional for the BA-225/225D/125D

Antistatic treatment


The glass panes of the breeze break are coated with transparent evaporated metal to block outside static electricity.



IR sensor

Daily/periodic balance check support (for the BA-T series)

The balance helps implement standard operating procedures (SOP) for quality and performance verifications without trouble. The daily check is to confirm that the balance is in good condition, whereas the periodic check is to regularly inspect the balance basic performance. No comprehensive knowledge or experience is required to implement these checks. Simply follow the displayed procedures step-by-step (or skip items if not needed for your laboratory).

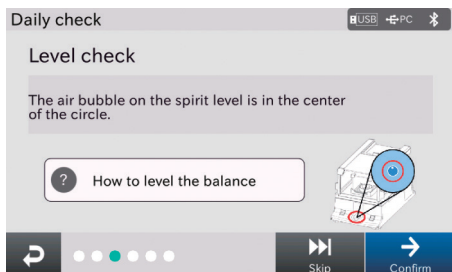


Daily check item: Level check

Other daily check items include external condition check, weighing pan check, and accuracy check.

Periodic check item: Eccentricity measurement


Other periodic check items include calibration test, sensitivity adjustment, repeatability measurement (with pre-loading), and sensitivity test.

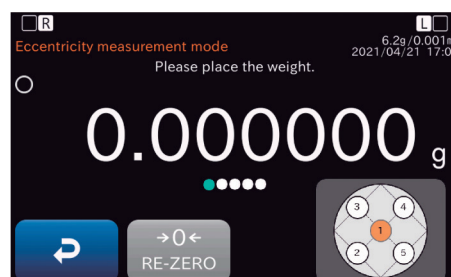

It is possible to set the balance to remind you of these checks as you prefer. Meanwhile, the results can be printed out or saved to a USB flash drive in a PDF report format for documentation and compliance.

Minimum weight setting and alert*2

You can set the minimum weight for the balance either by direct key input or performing a repeatability test, from which the balance automatically calculates*3 the minimum weight.

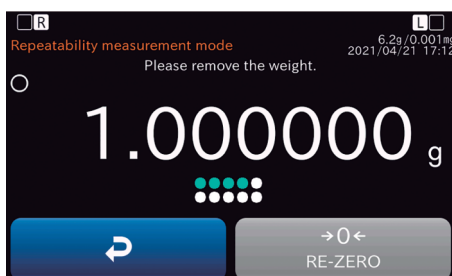

Repeatability measurement for minimum weight setting

The minimum weight is determined from the standard deviation of 10 repeated weighing results. The detailed calculation data can be output for your records.


*2 Unit of minimum weight setting is mg.

*3 By a tolerance of either 0.1% in accordance with the United States Pharmacopeia (USP), Chapter 41, or 1%, depending on your laboratory requirements

To ensure that the measured sample amount meets the minimum weight requirement, the balance can display an alert (message for the BA-T series) until the sample amount reaches the value entered as the minimum weight. You can also choose to set the balance to disable outputting of weighing data smaller than the minimum weight.

Minimum weight alert

You can select whether to include values near zero for comparison with the minimum weight.

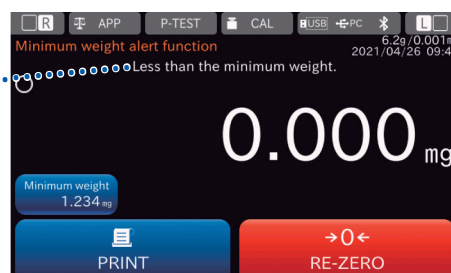

Advanced user access control (for the BA-T series*4)

For protection from unauthorized changes to balance settings/data, users can be classified into four levels (i.e. administrator, lab manager, supervisor, and operator), and the administrator can determine the extent of rights (i.e. change to settings, date/time setting, external sensitivity adjustment, and internal sensitivity adjustment) for each level.

User authorization				
	Change to settings	Date/time setting	Ext. sensitivity adjustment	Int. sensitivity adjustment
Administrator	Allowed	Allowed	Allowed	Not allowed
Lab manager	Allowed	Not allowed	Allowed	Not allowed
Supervisor	Allowed	Not allowed	Not allowed	Not allowed
Operator	Not allowed	Not allowed	Not allowed	Not allowed

User rights management

The administrator can register users as either lab managers or supervisors with usernames and passwords. Up to 100 users can be registered, including the administrator.*5

*4 For the BA series, the administrator can set passwords for up to 10 additional users. Similar (albeit simpler) user rights management is also possible using the function selection switches for these users as well as the administrator.

*5 Operators do not need a password.

History information (for the BA-T series)

The balance can display/save to a USB flash drive*6 log-in/log-out history, operation (changes to settings) history, and sensitivity adjustment history with date, time, username and other necessary information for later reference.

*6 The balance stores up to 1,000 data (which is then overwritten in order from the oldest data) for each history. It can only display the latest 100 data but save all data to a USB flash drive in CSV format.

Operation history			
	Date/time	User name	Item
1	2021/04/21 16:36:06	User 01	Buzzer sound
2	2021/04/21 16:36:13	User 01	Backlight Brightness
3	2021/04/21 16:36:16	User 01	Auto power OFF
4	2021/04/21 16:36:18	User 01	Auto power ON
5	2021/04/21 16:36:37	User 01	Auto door opening/closi
6	2021/04/21 16:37:07	User 01	Backlight Brightness

Operation history

GLP/GMP/GCP/ISO compliant output

Data necessary to comply with GLP/GMP/GCP/ISO can be added to sensitivity adjustment reports, calibration test reports, and series of weighing results.

Communication interfaces and connectors



- ① RS-232C (D-Sub 9P)
- ② Connector for the cable to the weighing unit
- ③ USB-A for USB flash drives*7
- ④ USB-B (mini-B) for connection with a PC
Using internal settings, you can toggle between Quick USB (HID) mode, which allows you to send weighing data to an application (e.g. spreadsheet) on a PC, and Virtual COM (CDC) mode for bi-directional communication. A USB cable (1.8 m) is supplied as standard.
- ⑤ Bluetooth®*7*8
- ⑥ Ethernet (TCP/IP)*7
- ⑦ External key input (jack socket) × 2
Separately sold foot switches can be connected to these sockets to operate the RE-ZERO, PRINT commands, or the automatic breeze break doors.
- ⑧ Connector for AC adapter

*7 For the BA-T series only

*8 The Bluetooth® function (GATT/HOGP) is currently enabled for the US, Canada and Japan only.

Specifications

	BA-6TE	BA-6DTE	BA-225TE	BA-225DTE	BA-125DTE
Capacity	6.2 g	2.1 g / 6.2 g ^{*i}	220 g	51 g / 220 g ^{*i}	51 g / 120 g ^{*i}
Readability	0.001 mg	0.001 mg / 0.01 mg	0.01 mg	0.01 mg / 0.1 mg	0.01 mg / 0.1 mg
Repeatability (standard deviation) ^{*ii}	0.0010 mg (for 1 g) 0.0030 mg (for 6 g)	0.0025 mg (for 1 g) 0.01 mg (for 6 g)	0.015 mg (for 50 g) 0.03 mg (for 200 g)	0.025 mg (for 50 g) 0.1 mg (for 200 g)	0.030 mg (for 50 g) 0.1 mg (for 100 g)
Minimum weight ^{*iii} (typical)	1.8 mg	4.0 mg	20 mg	32 mg	50 mg
Linearity	±0.010 mg	±0.010 mg / ±0.02 mg	±0.15 mg	±0.2 mg	±0.2 mg
Stabilization time (typical when set to FAST)	Approx. 10 secs	Approx. 10 secs / 10 secs	Approx. 7 secs	Approx. 7 secs / 3 secs	Approx. 7 secs / 3 secs
Internal weight ^{*iv}	Approx. 5 g		Approx. 100 g		
Sensitivity drift	±2 ppm/°C (10 to 30 °C/50 to 86 °F, when automatic self sensitivity adjustment is OFF)				
Operating environment	5 to 40 °C (41 to 104 °F), 85%RH or less (no condensation)				
Display unit	5-inch WVGA, TFT LCD color touch screen (resistive type) with two IR sensors				
Display refresh rate	5 times/sec or 10 times/sec				
Units of measure ^{*v}	mg (milligram), g (gram), oz (ounce), ozt (troy ounce), ct (metric carat), mom (momme), dwt (pennyweight), gr (grain), pcs (counting mode), and % (percent mode)				
Counting mode	Minimum unit mass	0.1 mg			
	Number of samples	10, 25, 50 or 100 pieces			
Percent mode	Minimum 100% reference mass	10.0 mg			
	% readability	0.01%, 0.1%, 1% (depends on the reference mass stored)			
Communication interface	RS-232C (D-Sub 9P), USB-A, USB-B (mini B), Ethernet (TCP/IP), External key input × 2, Bluetooth ^{®*vi}				
Applicable weights for calibration test/sensitivity adjustment	Any weight between 1 and 5 g		Any weight between 10 and 200 g		Any weight between 10 and 100 g
Weighing pan size	Ø25 mm		Ø85 mm		
External dimensions	Display unit	182 (W) × 138 (D) × 73 (H) mm			
	Weighing unit + breeze break unit	173 (W) × 305 (D) × 204 (H) mm		173 (W) × 305 (D) × 284 (H) mm	
	External ionizer	68 (W) × 129 (D) × 162 (H) mm			
Net weight	Display unit + weighing unit + breeze break unit	6.2 kg		6.7 kg	
	External ionizer	0.4 kg			
Power supply / consumption	AC adapter / approx. 36 VA				
Standard accessories	External ionizer × 1, Three sizes of aluminum analytical pans × 10 each (for the BA-6TE/6DTE), Tweezers for calibration weight × 1				

	BA-6E	BA-6DE	BA-225	BA-225D	BA-125D
Capacity	6.2 g	2.1 g / 6.2 g ^{*i}	220 g	51 g / 220 g ^{*i}	51 g / 120 g ^{*i}
Readability	0.001 mg	0.001 mg / 0.01 mg	0.01 mg	0.01 mg / 0.1 mg	0.01 mg / 0.1 mg
Repeatability (standard deviation) ^{*ii}	0.0010 mg (for 1 g) 0.0030 mg (for 6 g)	0.0025 mg (for 1 g) 0.01 mg (for 6 g)	0.015 mg (for 50 g) 0.03 mg (for 200 g)	0.025 mg (for 50 g) 0.1 mg (for 200 g)	0.030 mg (for 50 g) 0.1 mg (for 100 g)
Minimum weight ^{*iii} (typical)	1.8 mg	4.0 mg	20 mg	32 mg	50 mg
Linearity	±0.010 mg	±0.010 mg / ±0.02 mg	±0.15 mg	±0.2 mg	±0.2 mg
Stabilization time (typical when set to FAST)	Approx. 10 secs	Approx. 10 secs / 10 secs	Approx. 7 secs	Approx. 7 secs / 3 secs	Approx. 7 secs / 3 secs
Internal weight ^{*iv}	Approx. 5 g		Approx. 100 g		
Sensitivity drift	±2 ppm/°C (10 to 30 °C/50 to 86 °F, when automatic self sensitivity adjustment is OFF)				
Operating environment	5 to 40 °C (41 to 104 °F), 85%RH or less (no condensation)				
Display unit	Reverse backlit LCD (main characters: 11 segments, 17.8 mm height)				
Display refresh rate	5 times/sec or 10 times/sec				
Units of measure ^{*v}	mg (milligram), g (gram), oz (ounce), ozt (troy ounce), ct (metric carat), mom (momme), dwt (pennyweight), gr (grain), pcs (counting mode), and % (percent mode)				
Counting mode	Minimum unit mass	0.1 mg			
	Number of samples	10, 25, 50 or 100 pieces			
Percent mode	Minimum 100% reference mass	10.0 mg			
	% readability	0.01%, 0.1%, 1% (depends on the reference mass stored)			
Communication interface	RS-232C (D-Sub 9P), USB-B (mini B), External key input × 2				
Applicable weights for calibration test/sensitivity adjustment	5 g, 2 g, 1 g		200 g, 100 g, 50 g, 20 g, 10 g		100 g, 50 g, 20 g, 10 g
Weighing pan size	Ø25 mm		Ø85 mm		
External dimensions	Display unit	182 (W) × 138 (D) × 73 (H) mm			
	Weighing unit + breeze break unit	173 (W) × 305 (D) × 204 (H) mm		173 (W) × 305 (D) × 284 (H) mm	
	External ionizer	68 (W) × 129 (D) × 162 (H) mm		N/A	
Net weight	Display unit + weighing unit + breeze break unit	6.2 kg		6.7 kg	
	External ionizer	0.4 kg		N/A	
Power supply / consumption	AC adapter / approx. 36 VA				
Standard accessories	External ionizer × 1 (for the BA-6E/6DE), Three sizes of aluminum analytical pans × 10 each (for the BA-6E/6DE), Tweezers for calibration weight × 1				

*i Smart range function: Automatically switches between the precision and standard ranges. Changes back to full precision range when the RE-ZERO (tare) operation is implemented.

*ii Repeatability can worsen depending on the environmental conditions and operator skills.

*iii Pursuant to the United States Pharmacopeia (USP), Chapter 41


*iv The mass of the internal weight may change over time due to on-site environmental conditions and/or degradation with age.

*v Either tael (Singapore/HK jewelry/Taiwan) or tola can be added upon request.

*vi The Bluetooth[®] function (GATT/HOGP) is currently enabled for the US, Canada and Japan only.

Accessories

AD-1671	Anti-vibration table for balances
AD-1672/AD-1672A	Tabletop breeze break (large)
AD-1684A	Electrostatic field meter
AD-1687	Weighing environment logger
AD-1688	Weighing data logger
AD-8127	Compact printer
AX-BAT-31	Display cover for the BA-T series (5 pcs)
AX-BA-31	Display cover for the BA series (5 pcs)
AX-ION-25	External ionizer Also provided as standard for the BA-T series and the 0.001 mg models of the BA series
AX-BM-NEEDLESET	Discharge electrode units for the ionizer (a set of 4 pcs)
AX-IR-SWITCH	External IR switch Can be added to the ionizer if you prefer not to put a hand or sample close to the IR sensor of the ionizer.
AX-SW137-PRINT	Foot switch for PRINT (with connector)
AX-SW137-REZERO	Foot switch for RE-ZERO (with connector)

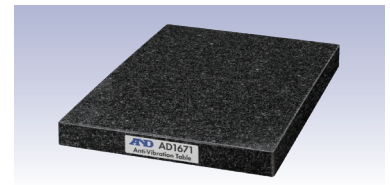

AD-1671
460 (W) × 400 (D) × 71 (H) mm




AX-ION-25 with AX-IR-SWITCH



Foot switches

For both series

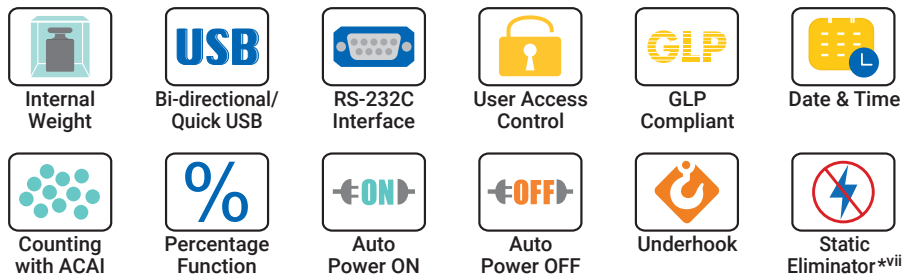

For the BA-T series only



For the BA series only



*^{vi} The Bluetooth® function (GATT/HOGP) is currently enabled for the US, Canada and Japan only.

*^{vii} Optional for the BA-225/225D/125D



Discover Precision

A&D Company, Ltd. (JAPAN)
URL: aand.jp

A&D Engineering, Inc. (USA)
URL: andonline.com

A&D Australasia Pty Ltd. (Australia)
URL: andaustralasia.com.au

A&D Instruments Ltd. (United Kingdom)
URL: andprecision.com

<German Sales Office>
URL: andprecision.com

A&D Korea Ltd. (South Korea)
URL: andk.co.kr

A&D Rus Co., Ltd. (Russia)
URL: and-rus.ru

A&D Instruments India (P) Ltd. (India)
URL: aandindia.in

A&D Sciencetech Taiwan Ltd. (Taiwan)
URL: aandd.com.tw

A&D Instruments Thailand Ltd. (Thailand)
URL: thai.andprecision.com