



## Increase Weighing Accuracy by Eliminating Electrostatic Charges

Electrostatic charges on weighing containers and samples can result in false readings or instability as they create attractive and repulsive forces. The ION-100A is designed to be used in conjunction with an OHAUS balance to neutralize electrostatic charges that can lead to inaccuracies.

#### Standard Features Include:

- Increases the Accuracy of Lab Measurements

  The ionizer neutralizes the electrostatic charge between the test tube and the balance, increasing the accuracy of measurements—particularly with low-volume samples.
- Fan-Free Technology Neutralizes Static with No Sample Disruption ION-100A's electrodes generate positive and negative ions that are delivered through fan-free technology that doesn't disturb fine samples.
- Compact and Versatile Design

  The ionizer can be adjusted for optimal alignment with the balance, and its small footprint takes up minimal space. The durable electrodes can be easily replaced and have an operating life of 15,000 hours.

## ION-100A Ionizer

## How does static affect your weighing results?


When the ambient relative humidity (RH) is below 45%, insulators such as plastics are typically charged with static electricity. This can cause errors of several milligrams, and impact weighing results. For example, powdery substances with a static charge might stick to to the edge of the test tube—making it difficult to weigh and to clean up.

#### **DC Corona Method**

Ionizer ION-100A uses the DC Corona Method to eliminate static charges. Two polarity electrodes continuously generate positive and negative ions. These ions are attracted to ions of the opposite polarity, neutralizing the sample within a charged environment.

Eventually, 99% of the static charge will be removed from the container. The most effective area is 10–40 cm away from the charged body.



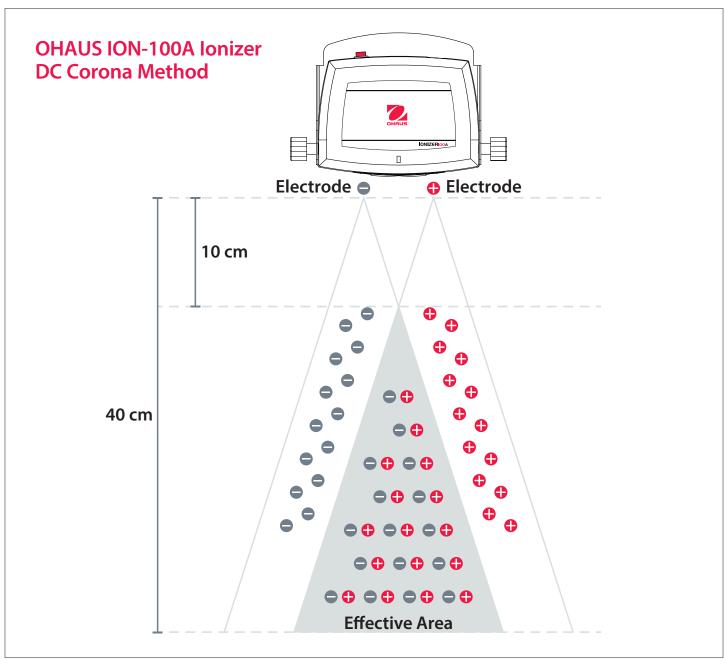


#### **DC Corona Method**


- Two electrodes work simultaneously to generate 1\*10<sup>6</sup>/cm<sup>3</sup> bipolar ions
- Covers a wide range of effective area lonization occurs in areas up to 40 cm away from the charged object
- Fan-free technology
   No need to worry about fine samples such as powders being blown away

#### **Conventional Method**

- One discharge needle generates a single type of positive or negative ions at intervals
- Static is removed most effectively in areas closest to the electrode only
- Requires a ventilator or an electronic fan to blow ions on to the sample

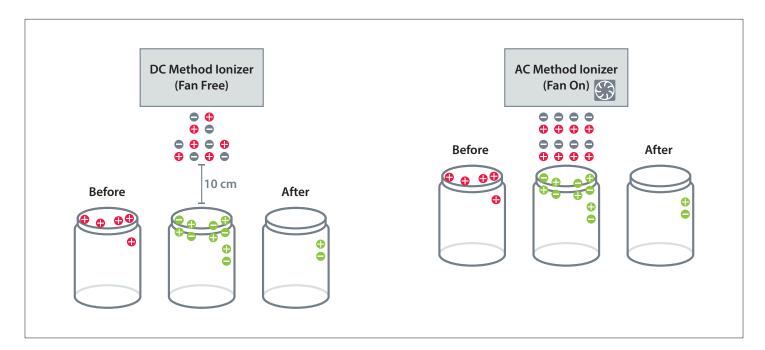

### **Dimensions**

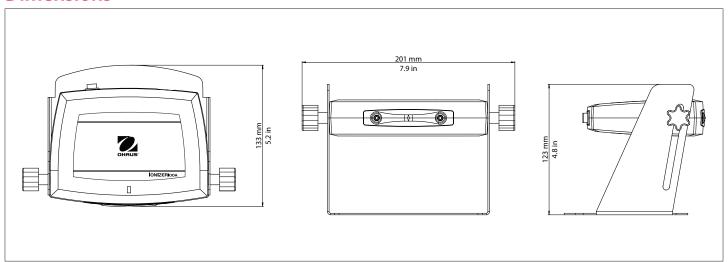

# ION-100A Ionizer

Model	ION-100A
Item Number	30130302
Power	12V == 0.5A
Ion Generation Method	DC Corona Discharge
Ionic Concentration	>1*10 <sup>6</sup> EA/cm <sup>3</sup> (10cm distance)
Effective Static Removal Range	100 to 400 mm from the outlet
Static Elimination Time (approx.)	1s/10 cm, 5s/20 cm (from ±1000V to ±100V)
Ozone Concentration	<0.05 ppm
Electrode Probes	Tungsten (φ2.5 mm/Length: 17.8 mm)(Durability: 15,000 hours)
Fan Speed	Fan-free
Ambient Conditions	Indoor use only
Altitude	Up to 2000 meters
Specified Temperature Range	0 °C to 40 °C
Humidity	Maximum relative humidity 80% for temperatures up to 31 °C (87.8 °F), decreasing linearly to 50% relative humidity at 40 °C (104 °F)
Power Supply Voltage Fluctuations	Up to ±10% of the nominal voltage
Installation Category	II
Pollution Degree	2
Operating Temperature Range	5 °C to 40°C / 41 °F to 104 °F
Net Weight	0.5 kg / 1.1 lb
Shipping Dimensions (W $\times$ H $\times$ L)	28 × 21 × 15 cm / 11 × 8.27 × 5.9 in
Shipping Weight	1.2 kg / 2.6 lb

### **Compliance**

- Product Safety: IEC 61010-1:2010, EN 61010-1:2010
- Electromagnetic Compatibility: IEC 61326-1:2012 (Class A emissions, General immunity), EN 61326-1:2013 (Class A emissions, General immunity), FCC Part 15 (Class A), Industry Canada ICES-001 (Class A)
- Environment: WEEE, RoHS

#### Accessory

OHAUS Europe GmbH Im Langacher 44

Im Langacher 44 8606 Greifensee Switzerland

e-mail: ssc@ohaus.com Tel: 0041 22 567 53 19 e-mail: tsc@ohaus.com Tel: 0041 22 567 53 20

www.ohaus.com

OHAUS Corporation is an ISO 9001:2008 manufacturer

ISO 9001:2008 Registered Quality Management System

CH16B001

